



Aerosart

Airfilter Cartridge for Industrial Applications

Membrane Filter Cartridges



Specifications

Materials

Filter Membrane:	PTFE
Support Fleece:	Polypropylene
Core:	Polypropylene
End Caps:	Polypropylene
O-Rings:	EPDM

Pore Size

0.2 µm

Available Sizes | Filtration Area

Size 1	10"	0.7 m ² 7.5 ft ²
Size 2	20"	1.5 m ² 16.1 ft ²
Size 3	30"	2.25 m ² 24.2 ft ²

Available Adapters

25, 27, 28

Packaging

6 cartridges per box

Operating Parameters

Max. Allowable Differential Pressure:	5 bar 72.5 psi at 20°C
	0.5 bar 7 psi at 134°C
Max. Allowable Back Pressure:	3 bar 43.5 psi at 20°C
	0.5 bar 7 psi at 134°C

Description

Aerosart high performance air filter cartridges can significantly reduce operating costs. The Aerosart is a high flow rate, low differential pressure, hydrophobic membrane filter. The unique single layer filter construction also reduces Blow-Down-time. Both the high flow rate and the short Blow-Down-time lowers the energy cost of air supply operations.

Applications

The Aerosart is designed for large-scale fermentation inlet and exhaust gas filtration.

Microbiological Safety

Aerosart filter cartridges have been tested and passed aerosol bacterial and viral challenge tests. Tests were conducted using MS-2 coli phages (NCIMB 10 108) and *B. subtilis* var. *niger* spores (NCTC 10073) at a challenge level of greater than 2.5×10^7 under worst case conditions of greater than 90% RH. No MS-2 coli phages or *B. subtilis* spores were detected on the downstream side of the Aerosart filter cartridges.

Performance

The unique single layer pleated filter construction of the highly hydrophobic PTFE membrane provides low differential pressures, excellent flow rates and the fastest blow down times of any gas service filter.

Long Service Life Time

The mechanical and thermal stresses experienced during steam in place sterilization pose the highest risk to any filter cartridge. In many cases, Aerosart filter cartridges will be used for more than the 120 steaming cycles. Tests have shown Aerosart cartridges to pass integrity tests with greater than 150 steaming cycles.

По вопросам продаж и поддержки обращайтесь:

Астана+7(7172)727-132, Волгоград(844)278-03-48, Воронеж(473)204-51-73, Екатеринбург(343)384-55-89,
Казань(843)206-01-48, Краснодар(861)203-40-90, Красноярск(391)204-63-61, Москва(495)268-04-70,
Нижний Новгород(831)429-08-12, Новосибирск(383)227-86-73, Ростов-на-Дону(863)308-18-15, Самара(846)206-03-16,
Санкт-Петербург(812)309-46-40, Саратов(845)249-38-78, Уфа(347)229-48-12

sst@nt-rt.ru || sartorius.nt-rt.ru

Specifications

Regulatory Compliance

Qualified for retention of aerosolized bacterial spores and viruses (coli-phages) in air

Non pyrogenic according to USP Bacterial Endotoxins

Passes USP Plastics Class VI Test

Non fiber releasing according to 21 CFR

Sterilization


In-Line Steam Sterilization

134°C, 20 min. at a maximum differential pressure of 0.5 bar | 7 psi

Sterilization Cycles

Minimum of 150 In-Line Sterilization cycles

Air flow rates Aerosart

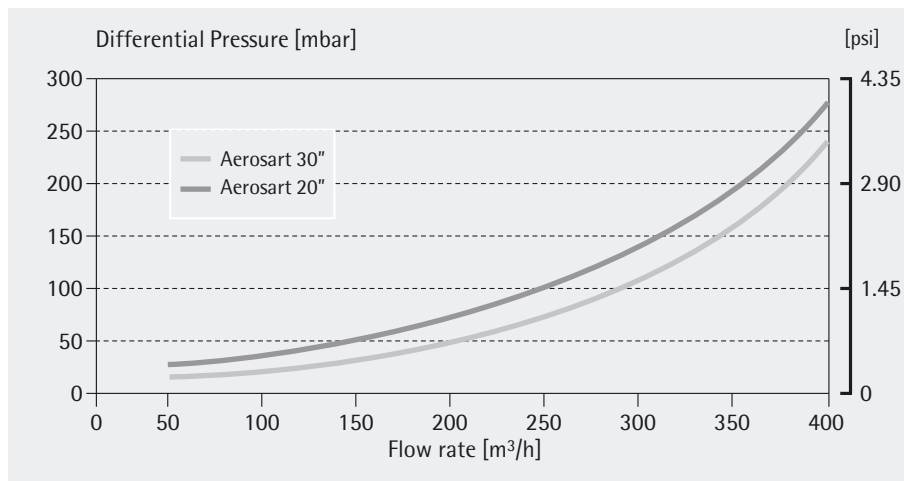

Air flow rate for Aerosart filter cartridges (0.2 µm) in relation to the filter cartridge heights at atmosphere pressure condition.

Order Information

Order Code	Pore Size [µm]	Size	Pack Size [Pieces]
5152507T1---E--C	0.2	10"	6
5152507T2---E--C	0.2	20"	6
5152507T3---E--C	0.2	30"	6
5152707T1---E--C	0.2	10"	6
5152707T2---E--C	0.2	20"	6
5152807T1---E--C	0.2	10"	6
5152807T2---E--C	0.2	20"	6

По вопросам продаж и поддержки обращайтесь:

Астана+7(7172)727-132, Волгоград(844)278-03-48, Воронеж(473)204-51-73, Екатеринбург(343)384-55-89,
Казань(843)206-01-48, Краснодар(861)203-40-90, Красноярск(391)204-63-61, Москва(495)268-04-70,
Нижний Новгород(831)429-08-12, Новосибирск(383)227-86-73, Ростов-на-Дону(863)308-18-15, Самара(846)206-03-16,
Санкт-Петербург(812)309-46-40, Саратов(845)249-38-78, Уфа(347)229-48-12

sst@nt-rt.ru || sartorius.nt-rt.ru